

## **PHYSITEMP INSTRUMENTS INC**

154 Huron Avenue

Clifton NJ 07013

Tel: 1-800-452-8510

Tel: 973-779-5577

Email: info@[physitemp.com](mailto:info@physitemp.com)

### **Instructions for Sterilization of T204E Flexible Microprobes**

#### **Cleaning before sterilization:**

Wipe with alcohol or cleaning solution

Note: The plug should never be immersed in liquid. It is sealed with epoxy to prevent moisture entering during normal use and wiping, but has not been tested for immersion.

**AUTOCLAVING:** Sterilize at: 134°C for 3 minutes at 2.2 bar  
or 121°C for 15 minutes  
or 115°C for 30 minutes